


JEOL 502 aka Ripley

How to open your file on a workstation


Last update 3/8/16

Ripley (spectrometer): 18.60.3.245


1. Click on the Delta icon. Click on the folder icon.

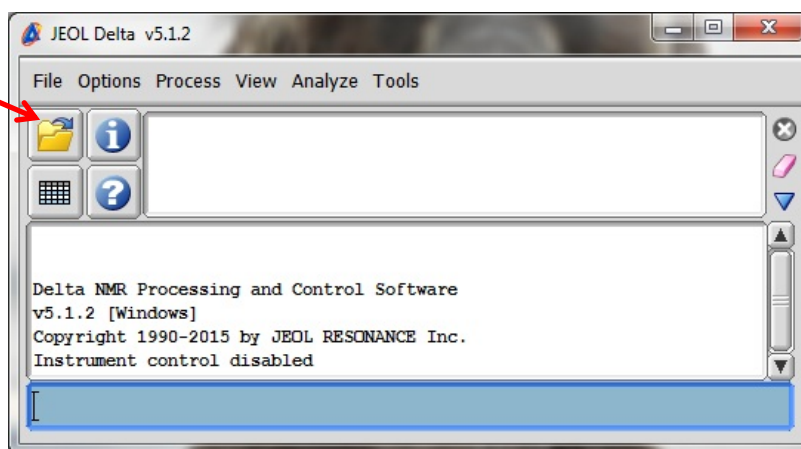


2. You will get this window. Any listed files are ones already saved locally. Click on Data Servers.

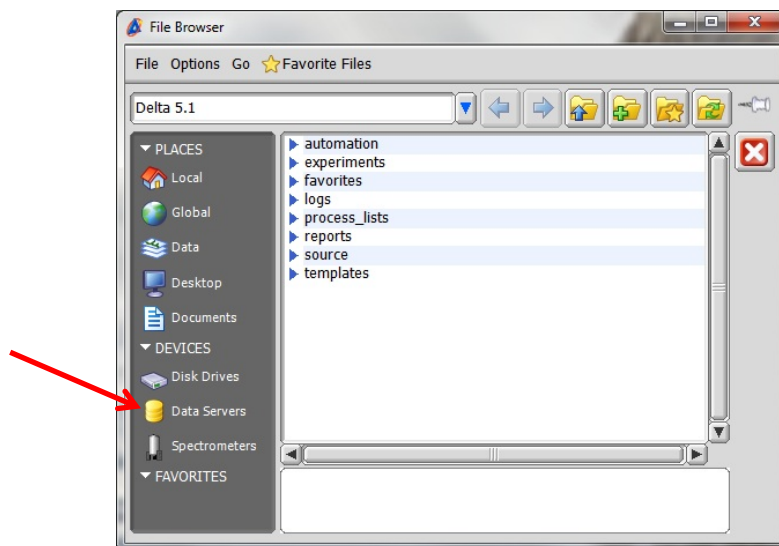


3. You will get this window. Select Ripley.

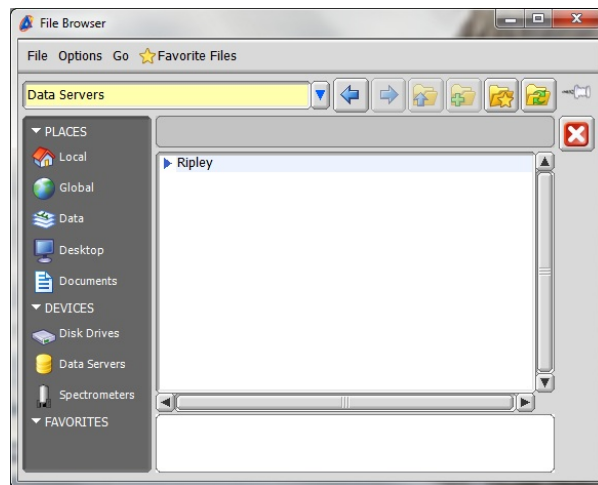


4. Click **Connect As...**

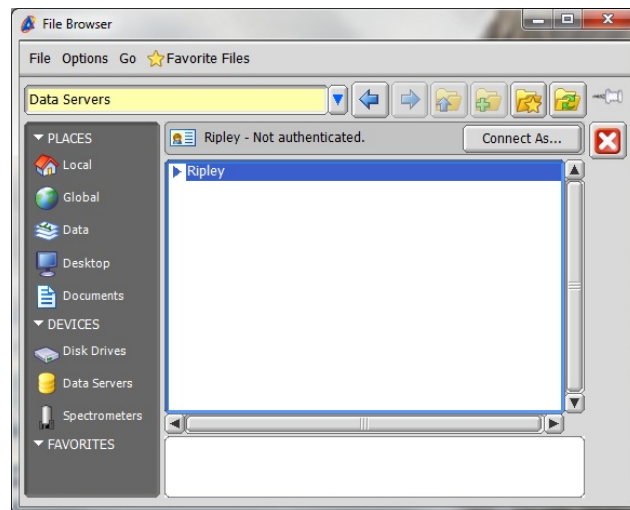


5. Log yourself in.

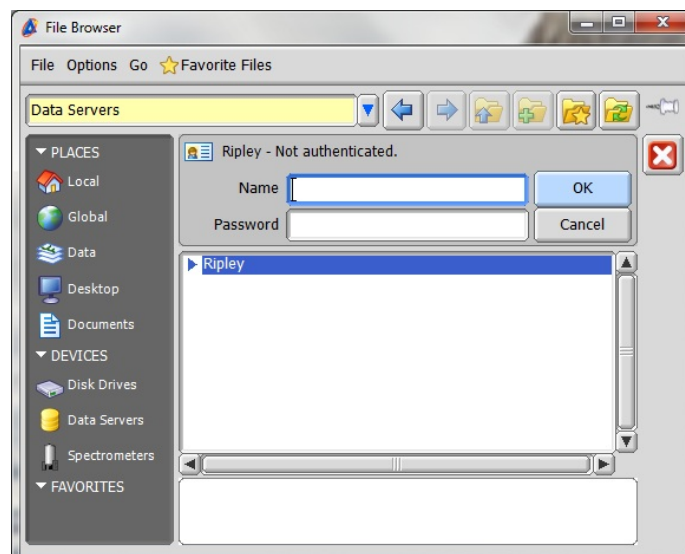


6. Next you get a list of your files on the spectrometer data server.

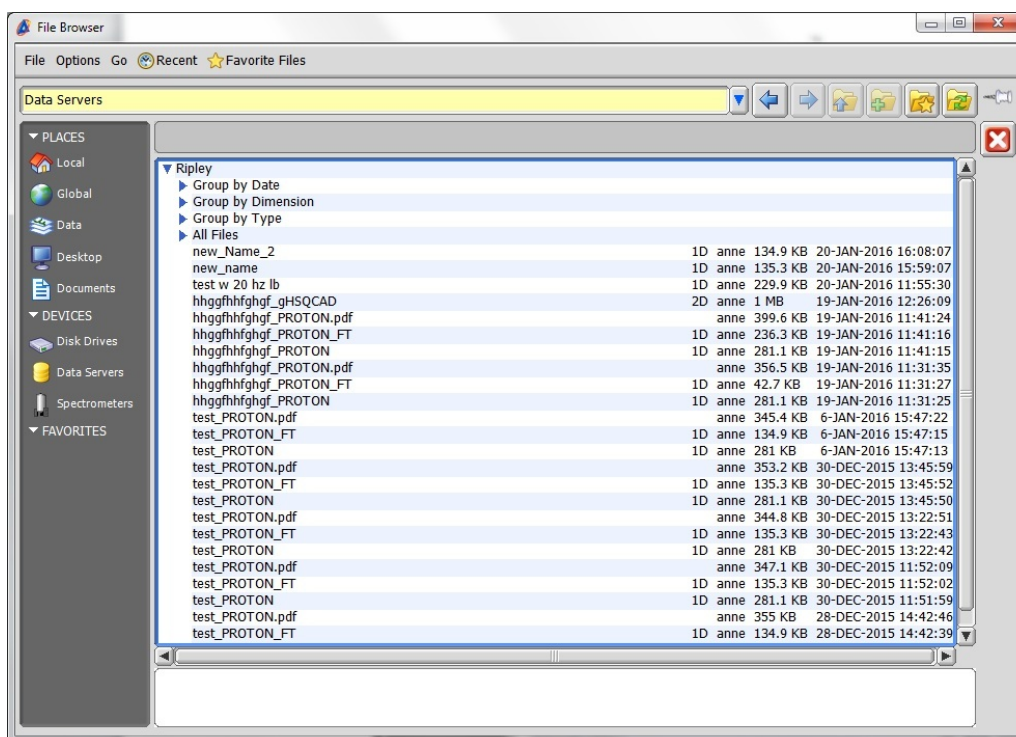


7. Double click on the file you want to work up (or click on the **green arrow**). Your file should open.

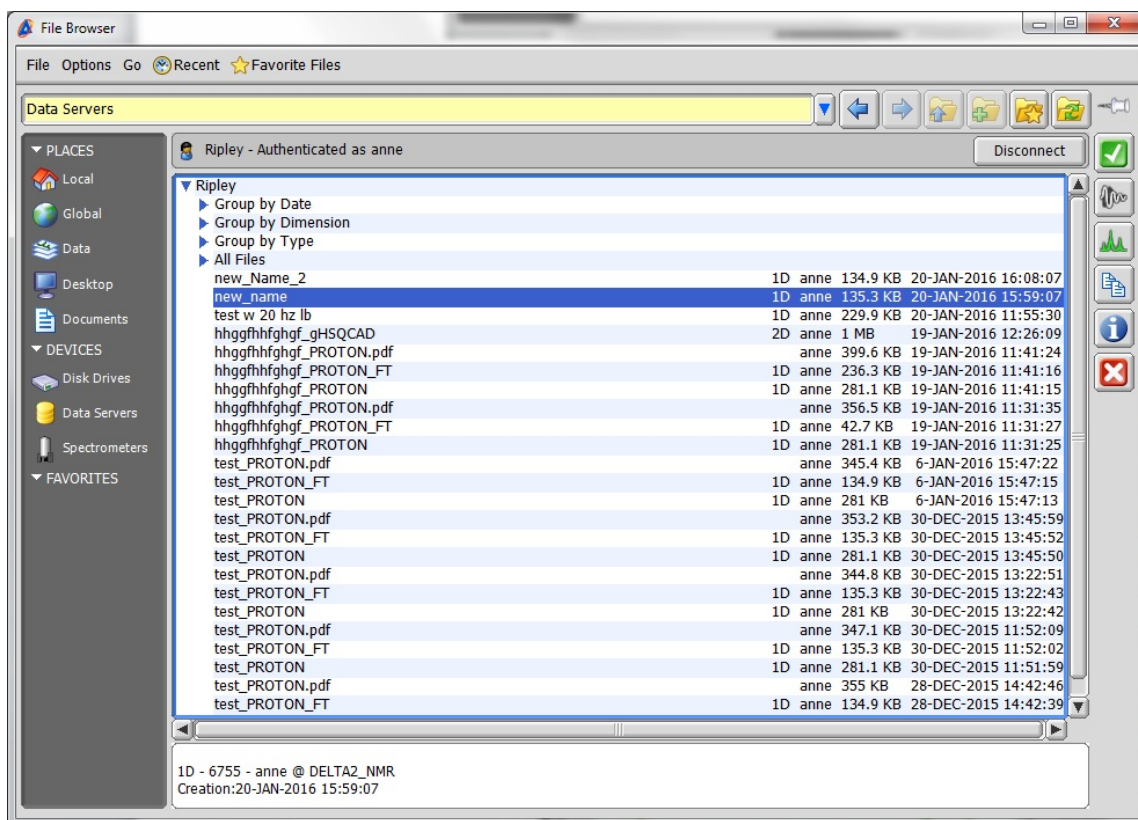

8. You can save your changes locally (on the workstation). Go to File → Save As...

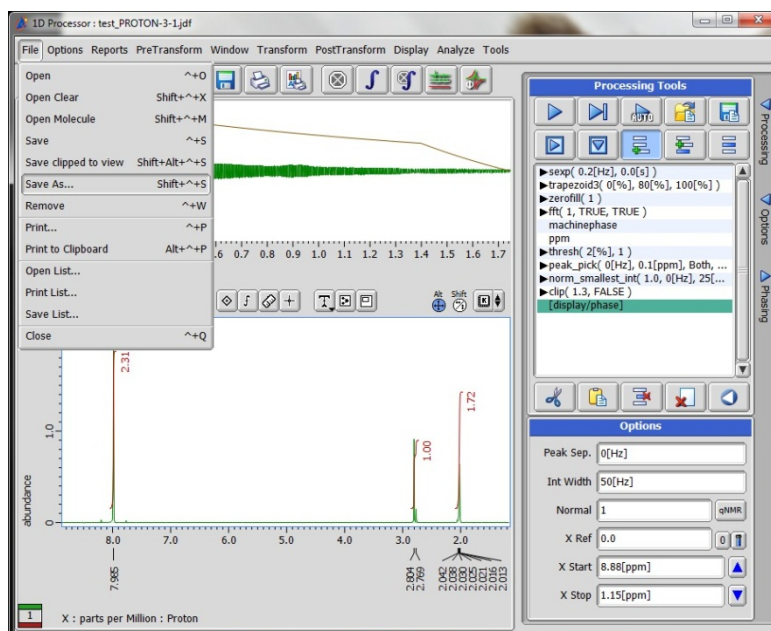


9. Rename it and your processed data will be saved locally. The raw file is still on the spectrometer data server, and can be accessed from another workstation.

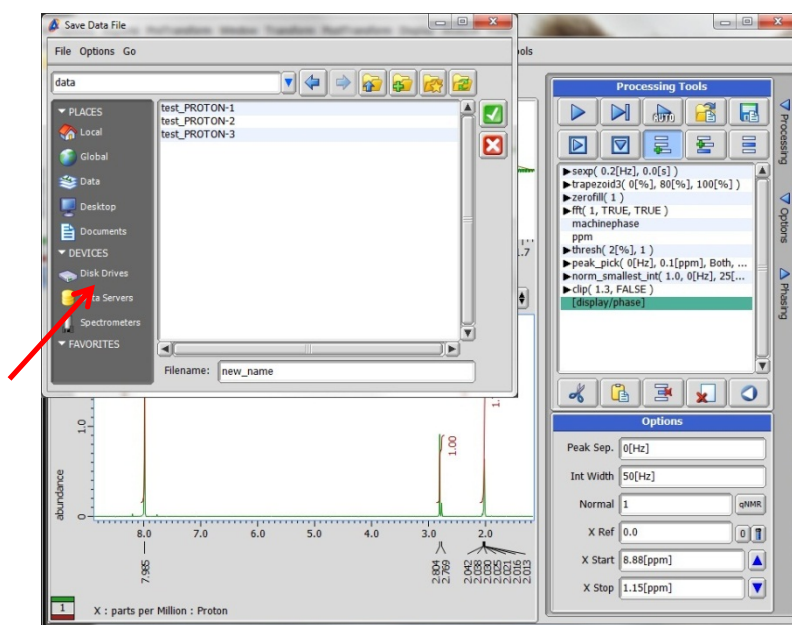


If you want to save a file to a USB stick or flash drive, select Disk Drives under DEVICES.


10. If you want to save your changes to the spectrometer's data server, rename the file in the **Save As...** window.

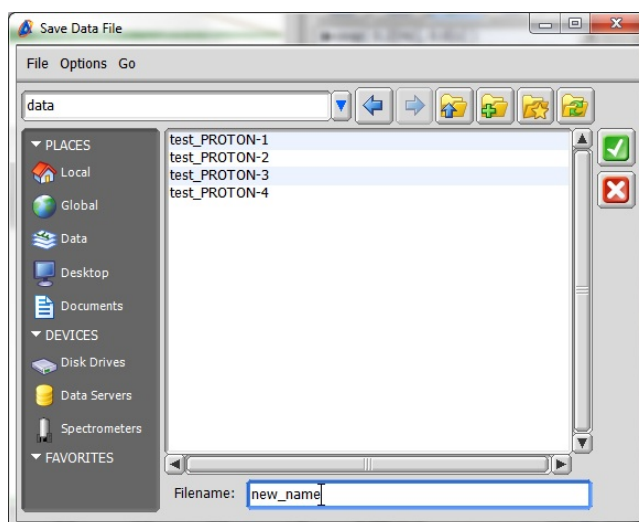


11. Go back to this window and click the folder again.

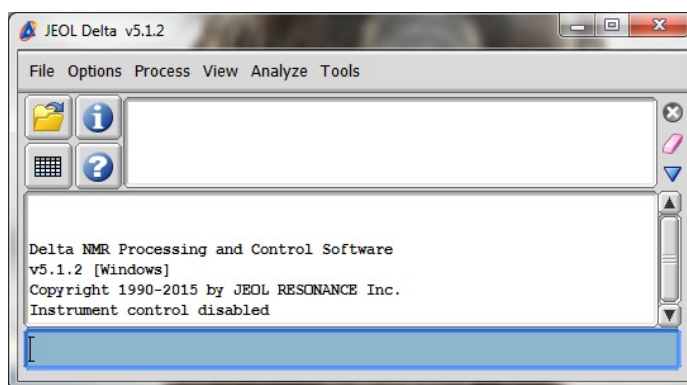



12. The newly named file will appear in Data. Select the file you want to save.

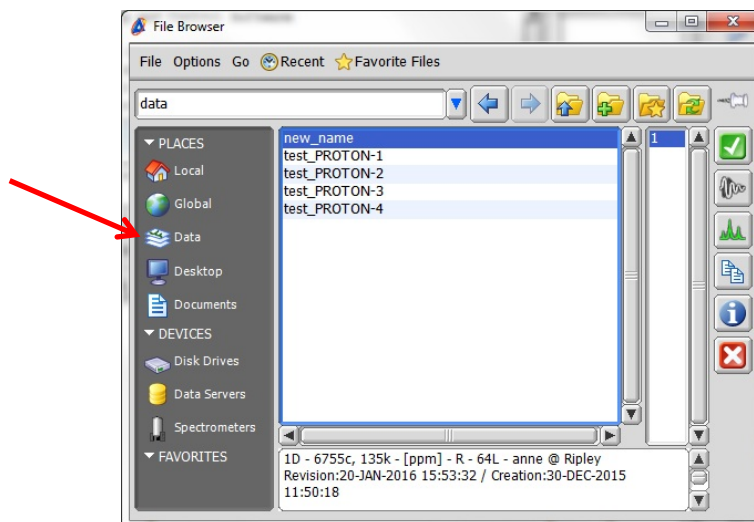

13. File → Upload File...

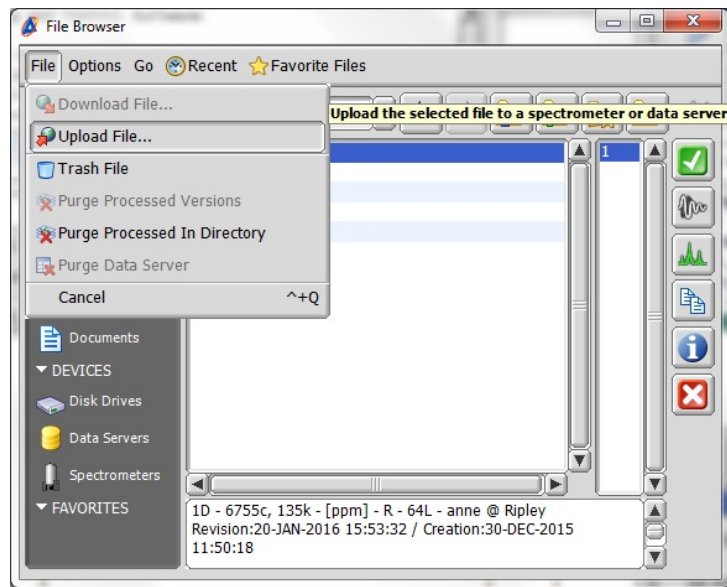

14. Select the Data Server, then Ripley, and click the **green arrow**.

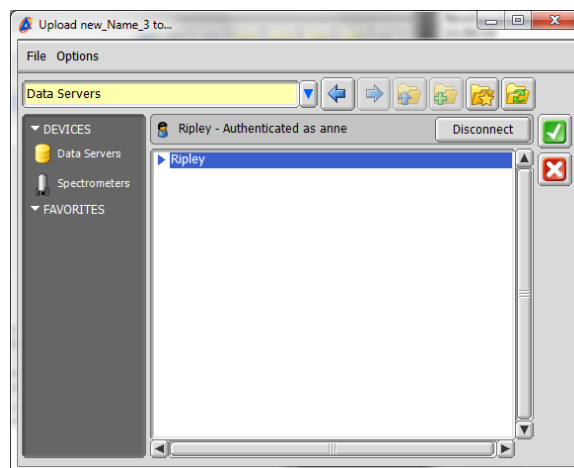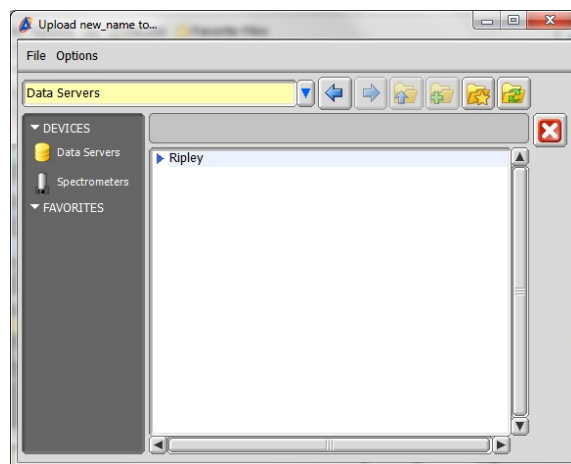

Your file should be saved back to the spectrometer's data server.

Disconnect when done, and close Delta.